16214

Directions for Use
Stainless Steel Syringe Filter Holder

1. Disassemble the syringe filter holder and carefully clean the parts before using for the first time and after each use.

2. After cleaning, rinse the parts; use distilled water for the last rinse.

3. Place the parts in a drying oven to dry or blow-dry with compressed air.

4. Position a membrane filter in the recess of the base using forceps.

5. If you need to use a prefilter or glass fiber filter, center it on the membrane filter.

6. Keep the filter holder horizontal when screwing the parts together.

7. Place the filter holder in the plastic locking tool and tighten the assembled halves.

8. Remove the syringe filter holder from the locking tool.

9. Attach a hypodermic needle to the male Luer slip of the syringe filter holder, and wrap the needle in aluminum foil.

10. Autoclave the prepared filter holder for 30 minutes at 121°C. Check the operating temperature of the autoclave from time to time with a maximum recording thermometer.

11. When the autoclaving time is up, allow pressure equilibration to take place in a non-programmed autoclave by cooling.

12. Allow the autoclaved syringe filter holder(s) to cool off and retighten with the locking tool.

13. If cellulose acetate membrane filters 111.. are used, the filter holder(s) can be sterilized with dry heat at 180°C for 2 hours instead of autoclaving.
14. Draw up the liquid to be filtered into the syringe.

15. Attach the sterile syringe filter holder to the Luer lock connector of the syringe.

16. Remove the aluminum foil from the needl before filtering.
- For the purpose of product development we reserve the right to make changes without notice.

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Specifications subject to change without notice. Printed and copyrighted by Sartorius Stedim Biotech GmbH
Sartorius Stedim Biotech GmbH